

Lithium Ore

DIGESTION OF LITHIUM ORE USING COLDBLOCK™ DIGESTION TECHNOLOGY

Author: Lorie-Anne Doig

Introduction

This application note will focus on the digestion of lithium ores using ColdBlock™ Digestion CB12L Technology.

Method

Triplicate samples were digested using the following method:

- 1.0g sample was weighed and placed into a ColdBlock™ Digestion vessel.
- 25mL Aqua Regia was added (3:1 HCl:HNO₃)
- Sample was digested at 100 % power for 10 minutes; chiller temperature was set to -5°C.
- Sample was allowed to cool then was normalized to 50mL using 2% HNO₃.

Instrument

ColdBlock™ Digestion CB12L Technology.

General

This procedure is specific for the sample digested and may need modification for different samples to achieve the desired result.